



# BD™ Sterile Pack Media Products

## PERSONNEL AND ENVIRONMENTAL MONITORING



### Environmental monitoring of your ISO Class 7 sterility suite:

- Upon entry take active air sample.
- Sanitize hood.
- Expose 2 settling plates (100 mm) on hood throughout testing.
- Take 1 RODAC™ plate - sample floor and wall per SOP.
- Take swab sample for irregular surfaces.
- Monitor personnel at exit.

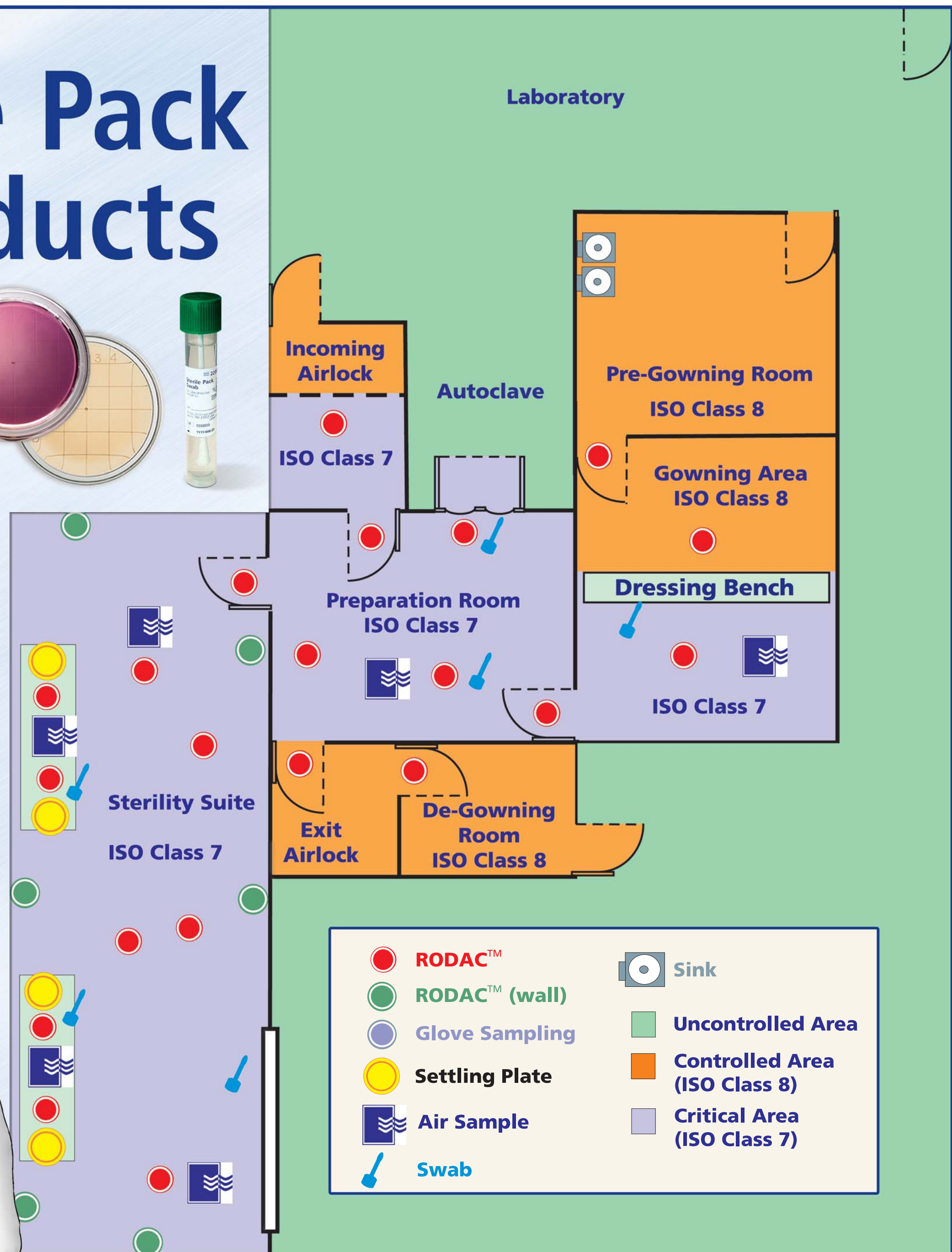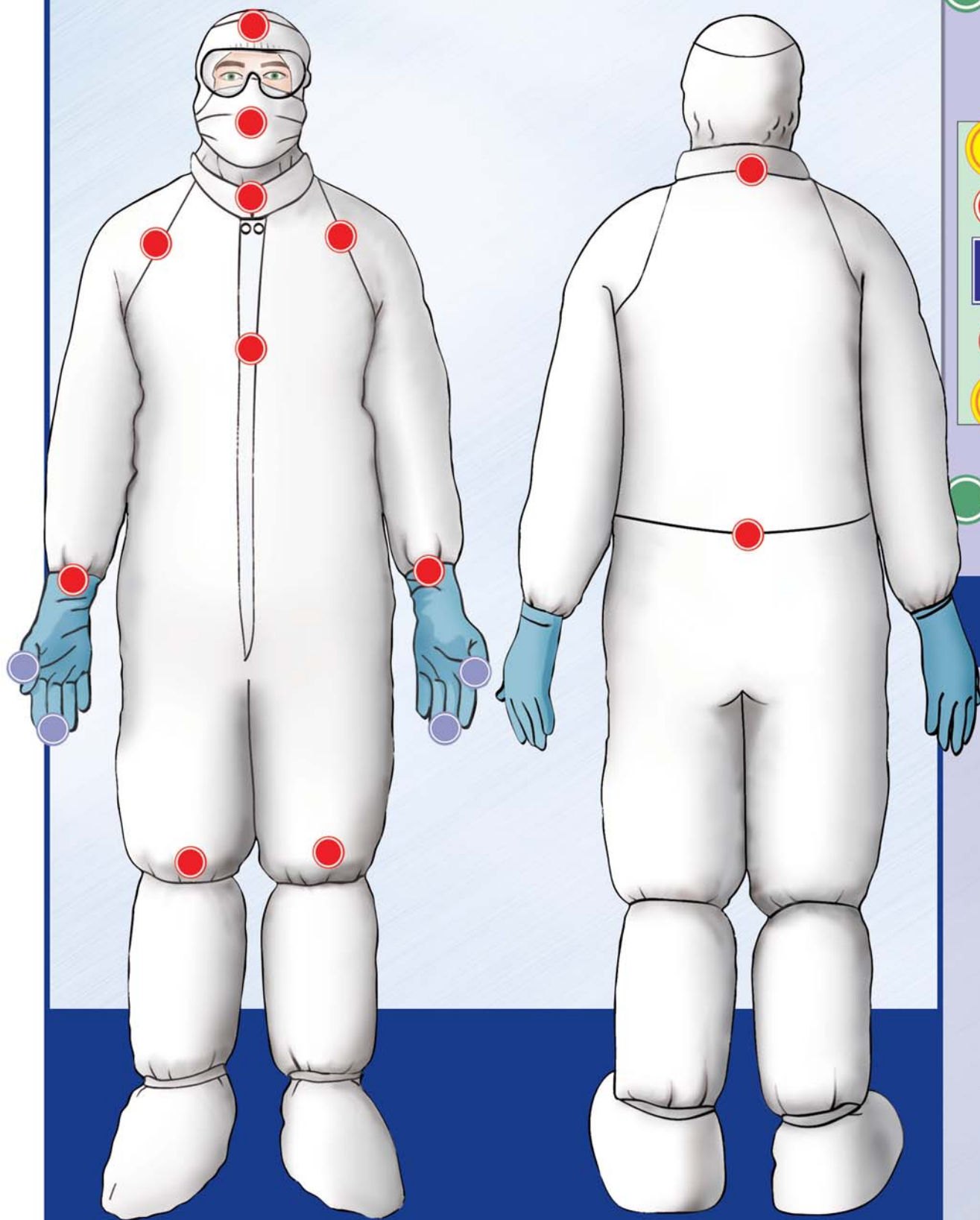

Description	Unit/Quantity	BD Cat #
<b>BD BBL™ Sterile Pack RODAC™ Plates for Personnel and Surface Sampling*</b>		
D/E Neutralizing Agar	Package of 10	221232
	Carton of 100	222209
Trypticase™ Soy Agar	Package of 10	221238
with Lecithin and Polysorbate 80	Carton of 100	222208
Sabouraud Dextrose Agar	Package of 10	221233
with Lecithin and Polysorbate 80	Carton of 100	292351
<b>BD BBL™ Finger Dab™ Plates for Personnel Monitoring*</b>		
Trypticase™ Soy Agar	100 mm Package of 10	292273
with Lecithin and Polysorbate 80	100 mm Carton of 100	292271
	150 mm Package of 5	292305
<b>BD BBL™ Sterile Pack Settling Plates for Active and Passive Air Sampling*</b>		
Trypticase™ Soy Agar	100 mm Package of 10	221236
	100 mm Carton of 100	222205
	150 mm Package of 5	221237
	150 mm Package of 45	222206
<b>BD Sterile Pack Swab for Surface Sampling</b>		
Sterile Rinse Solution, 10 mL	Pack of 200	220518



**BD Diagnostics**  
 7 Loveton Circle  
 Sparks, MD 21152-0999 USA  
 800.638.8663  
 www.bd.com/ds

\*Please see BD catalog for complete BD Sterile Pack Media product offering.

BD, BD Logo and all other trademarks are the property of Becton, Dickinson and Company. ©2004 BD. 1-2406 July 2004 Printed in the USA