

BD Sterile Pack Swabs

For Environmental Monitoring



Maximum sterility with double wrapped packaging, gamma-irradiation and performance-validation.



Helping all people
live healthy lives

BD Sterile Pack Swabs For Environmental Monitoring

Maximum sterility with double wrapped packaging, gamma-irradiation and performance-validation.



DON'T TAKE CHANCES



that your environmental samples could be contaminated by a less than sterile swab. Wipe out uncertainty with BD Sterile Pack Swabs – the first pre-filled swab/rinse solution set with a Sterility Assurance Level (SAL) of 10^{-6} , an industry standard! BD Sterile Pack Swabs are ideal for the stringent sterility requirements of surface sampling in clean rooms and isolators at pharmaceutical and medical device manufacturing facilities. And because these swabs are designed for hard-to-reach places, BD Sterile Pack Swabs easily fit into equipment recesses, nooks and crevices.

PRODUCT FEATURES

- Packaged 50 to a box, with 10 single wrapped swabs in each double wrapped pouch
- Polypropylene tube is pre-filled with a 10 mL rinse solution
- Simple sampling and transport—easy-to-use swab is fixed to the screw cap
- Complies with the new ISO guidelines
- Each lot comes with a Certificate of Analysis
- ATP-free, Dacron™ swab
- Saves time, no need to create your own media and swab/solution set in house
- Convenient and cost-effective solution
- One year shelf-life, at room temperature storage


PROCEDURAL STEPS



1. Open double wrapped exterior and peel open single wrapped pouch.



2. Twist cap and remove swab from tube.

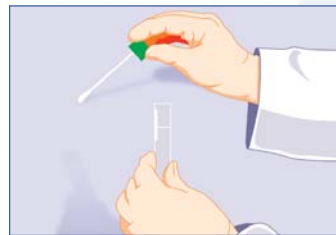

3. Swab sampling area.



4. Twist cap onto tube and transport to the laboratory for testing.

ORDERING INFORMATION

Cat. Number	Description	Qty/Pkg.
220518	BD Sterile Pack Swabs	each



BD Diagnostics

7 Loveton Circle
Sparks, MD 21152-0999 USA
800.638.8663
www.bd.com/ds

11 rue Aristide Bergès
38800 Le Pont de Claix, France
Tel: 33.4.7668.3636

2771 Bristol Circle
Oakville, Ontario
Canada L6H 6R5
Tel: 800.268.5430

30 Tuas Avenue 2
Singapore 639461
Tel: 65.6861.0633

Monte Pelvoux 111, 9th Floor
Col. Lomas de Chapultepec
11000 México D.F.
Tel: 52.55.59.99.82.00

Rua Alexandre Dumas 1976
04717-004 São Paulo, S.P. Brazil
Tel: 55.11.5185.9833

Akasaka DS Building
5-26 Akasaka 8-chome
Minato-ku
Tokyo, 107 Japan
Tel: 81.24.593.5405